


Main Menu and Navigation Flowcharts


AD Series

(Automatic Zero IZSM) Press T when Auto is displayed.

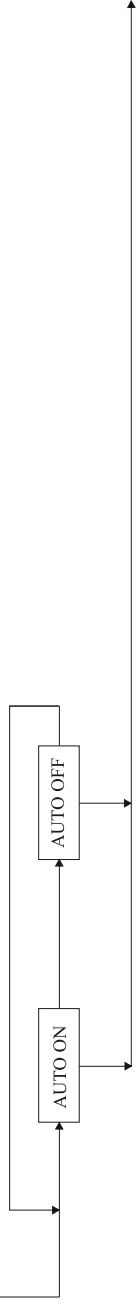


Press T when the desired selection is displayed.

(GLP) Press T when GLP is displayed.

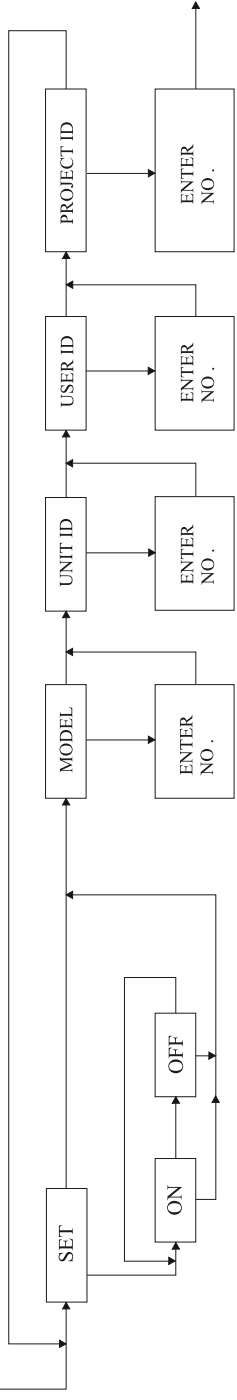


(Print) Press T when PRINT is displayed.

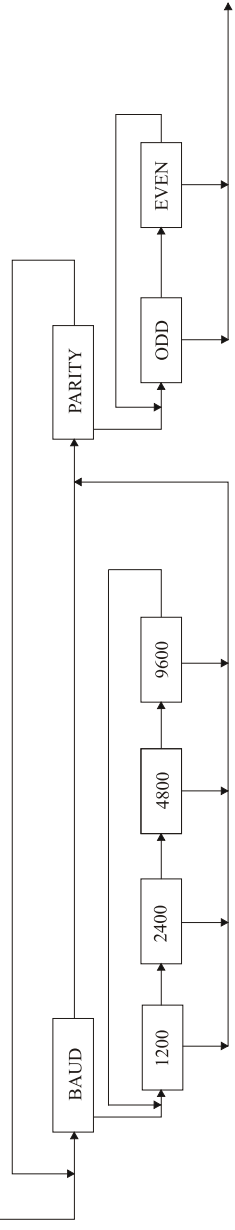

Press T when the desired selection is displayed.

(Time and Date) Press T when SET TD is displayed.

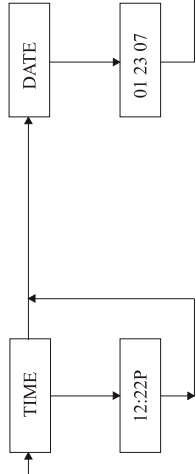

Press the →0< (NO) key to change a digit and the T (YES) key to accept your selection.


(Master Reset) Press T when RESET is displayed.



Press the T (YES) key to confirm the Master Reset or →0< to Cancel and Abort.


To enter menus, accept commands or confirm options use the “T (YES)” key. To exit, abort or cancel use the “C”.

(Main Menu) Press M key to enter the Main Menu

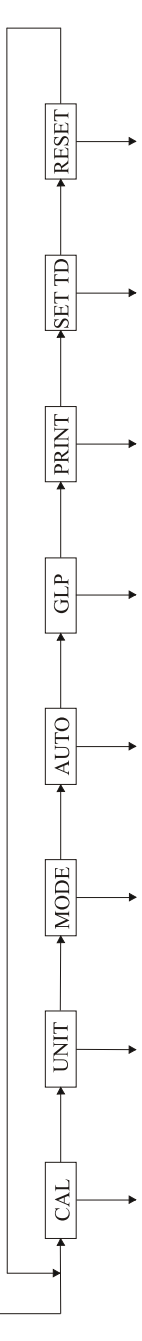

Press T when desired submenu is displayed.

(CALIBRATION) Press T when CAL is displayed.

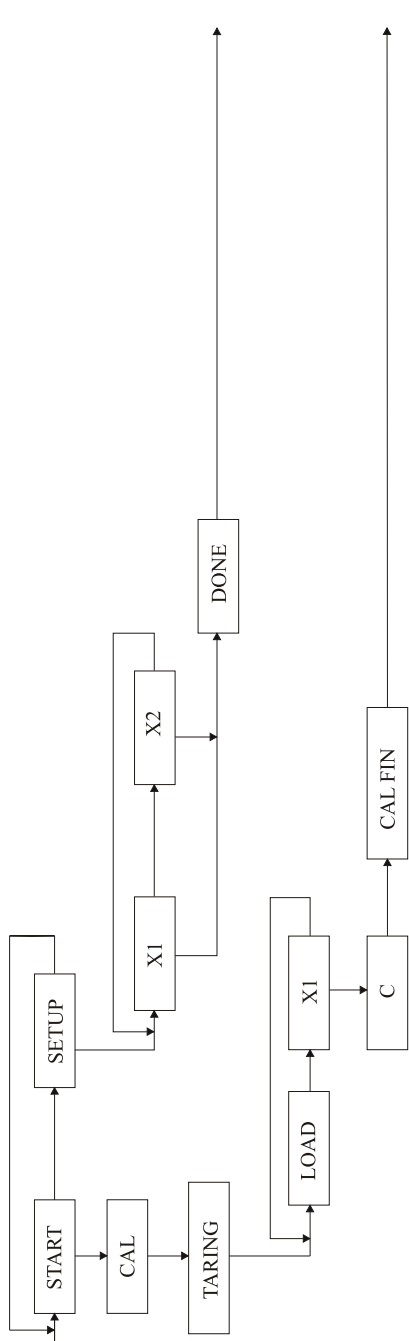

(UNIT) Press T when Unit is displayed

